

DigiGenius Gel Image Capture-Quick Guide

1 Turn on

Select



- ➔ Turn on the camera (make sure the battery is charged)
- ➔ **N.B.** If the Canon Powershot software launches automatically please complete the following step before using the software
- ➔ Click on Start/All Programmes/Canon utilities/camera window/remote capture DC3/remote capture DC3

2 Capturing an image


Select



- ➔ From the Canon camera window select Remote shooting tab
- ➔ Press start
- ➔ Select directory you wish to store the images in

3 Turn on transilluminator

Select

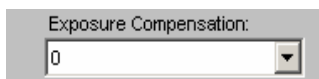

- ➔ If bands on gel are faint then increase the exposure compensation from the shooting settings menu
- ➔ The image is now live

4 Zoom setting

Select



- ➔ To adjust the level of zoom use the slider bar to zoom the camera in or out

5 Grab image

Press



- ➔ Then turn off the camera

6 Save and export

- ➔ Export as TIFF or BMP
- ➔ Images are saved as JPEG files

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
DGIC0210