

# GelVue Dual UV Transilluminator

## Description

GelVue Dual UV Transilluminator is an ultra violet light box with two UV wavelengths, for use in gel imaging.


## Model

Model	Filter size	Wavelength (nm)	
GVLM20	20 x 20	365 and 302	L = Long = 365
GVLM30	20 x 30	365 and 302	
GVLM25/30	25 x 30	365 and 302	M = Medium = 302
GVLS20	20 x 20	365 and 254	S = Short = 254
GVLS30	20 x 30	365 and 254	
GVLS25/30	25 x 30	365 and 254	
GVSM20	20 x 20	254 and 302	
GVSM30	20 x 30	254 and 302	
GVSM25/30	25 x 30	254 and 302	

**Warning: UV radiation can be dangerous for unprotected eyes and skin, therefore, appropriate precautions must be taken. UV protective goggles or a face shield must be worn. Please contact our sales office for further details and accessories.**

## Operating Instructions

- 1) Check that mains cable is correctly earthed and that plug is fitted with 5 amp fuse.
- 2) Check voltage supply requirements on the side of the unit.
- 3) Check that there is nothing loose inside the unit.
- 4) Note the switch arrangement is as shown in the diagram below. Select the wavelength switch of choice. Make sure that the On/Off switch is in the Off position.

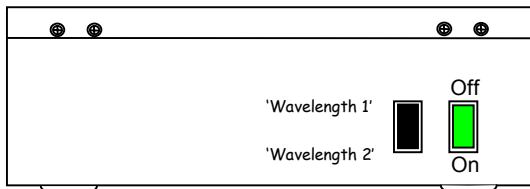

'Wavelength 1' = L/L/S if

'Wavelength 2' = M/S/M

- 5) Plug the mains input to the unit and supply. **Do not switch the unit on yet.**
- 6) Make sure that the area is safe, ie, safety precautions must be observed.
- 7) If safe, switch the unit on and ensure the switch lamp illuminates and the lamps in the unit illuminate evenly. Note only half the lamps should illuminate (every other one).
- 8) It is our recommendation to switch the unit off if not in use.

## If the unit fails to light up

Check that the mains supply is present. Check that the On/Off switch is switched to On.

If unit still fails to light, check the fuse in the unit and plug.

If no fault can be found, contact our service department for advice and assistance.



### CONTACT SYNGENE:

EUROPE  
BEACON HOUSE  
NUFFIELD ROAD  
CAMBRIDGE  
CB4 1TF  
Tel: +44 (0)1223 727123  
Fax: +44 (0)1223 727101  
Email: sales@syngene.com  
Email: support@syngene.com

[www.syngene.com](http://www.syngene.com)

IN USA:  
5108 PEGASUS COURT, SUITE M  
FREDERICK  
MD 21704

Tel: 800 686 4407/301 662 2863  
Fax: 301 631 3977  
Email: ussales@syngene.com  
Email: support@syngene.com