

# GeneGnome HR Image Capture-Quick Guide

## 1 Turn on

➔ Press the black button at the rear of the instrument

## 2 Start GeneSnap software

➔ Each user can enter a name or title for their own individual user preferences

## 3 Open drawer

Select



Open



Closed

➔ Place your sample in between the scored 8 position arrows present on the edges of the drawer frame. Then gently push the drawer to close

## 4 Choose appropriate settings

Select



Press



Select




➔ Select the classic tab  
➔ When the door is shut the button will turn blue and the image is no longer live

➔ Press the blue button to start

➔ Select series capture to either capture a single frame or a series of images

➔ Using the drop down menu select the number of images you wish to capture in the series

➔ Set exposure time and click ok

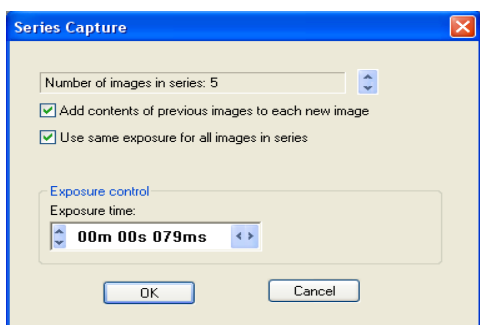

5

## Save

Click



- ➔ Choose from Save As
- Or
- ➔ File/Export image to TIF, BMP, or JPEG

6

## Print

Click



- ➔ Sends image to printer

7

## Help menu

- ➔ Select the help menu from the toolbar
- ➔ Use the drop down menu to pick a help topic

CONTACT SYNGENE:  
EUROPE:  
Tel: +44 (0)1223 727123  
Email: support@syngene.com

USA:  
Tel: 800 686 4407/301 662 2863  
Email: support@syngene.com

[www.syngene.com](http://www.syngene.com)  
GGHRClassIC0210