

GeneGnome HR Installation - Quick Guide

■ System components

1. GeneGnome HR (red mini darkroom)
2. Integral PC
3. Flatscreen monitor with keyboard and mouse

■ Installation

Connecting the components

1. Unscrew the cover plate on the left-hand side of the GeneGnome HR
2. Remove the cover plate to reveal the connectors beneath
3. Plug in the flat panel monitor, keyboard and mouse leads (the sockets are labelled with appropriate icons)
4. If a thermal printer has been supplied connect the printer cable to the appropriate socket
5. Refit the cover panel, slotting the cables through the cut-outs

CAUTION: Connect electrical power leads to each component but DO NOT SWITCH ON AT THIS TIME

■ Switching on

1. Ensure each mains lead is now connected to a live supply
2. Switch on the GeneGnome HR using the On/Off switch at the rear of the unit
3. Switch on the GeneGnome HR processor (hold the round button on the rear of the GeneGnome HR for 1-2 seconds). Also switch on the computer monitor and thermal printer (if connected)
4. The software is pre-loaded. You will need to accept the Windows licence

When the Windows operating system is fully loaded the GeneGnome HR system is ready to use

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
GGHRINST1209