

U:Genius Image Capture-Quick Guide

1 Turn on

- ➔ Press the round black button at the top of the back of the unit
- ➔ Position sample
- ➔ To turn on UV close door and check UV is on, on front of unit

2 Capturing an image

Click




- ➔ A live image will be displayed on the screen

3 Camera settings

Aperture ➔

Zoom ➔

Focus ➔

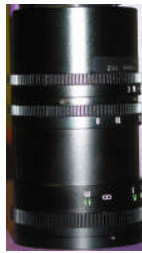

- ➔ Adjust Aperture, Zoom and Focus settings until a suitable image is displayed, preferably with sample filling the screen for maximum resolution
- ➔ N.B. Opening the aperture too far may result in areas of the sample being saturated generating a very bright image

4 Exposure settings

Manual settings

Click



Auto-exposure setting



- ➔ Click (+) to increase and (-) to decrease the exposure time
- ➔ Auto-exposure function sets the exposure time to a level that ensures no saturation of the image occurs. If you use this button image is captured and frozen

5 Grab image

Press (manual exposure only)



- ➔ Image will now be frozen

6 Image contrast slider bar



- ➔ The image contrast slider bar enables the appearance of the image to be changed
- ➔ Move the green/blue line to the left to lighten the image and move to the right to darken the image

7

Changing image settings

Zoom setting

Press



➔ Zoom controls

Use the (+) to zoom in and the (-) to zoom out

➔ When zoomed in you can 'pan' the image by dragging with your finger

Press



then press



Rotate setting

➔ Rotate

Rotate the image by touching a point on the image and move your finger on the screen

Press



then

Sharpen Smooth Invert



➔ Image processing

You can sharpen, smooth and invert image

Press



then



Annotate setting

➔ Annotate

To add and edit text annotations press the area of the image you want to annotate. To select text you can highlight the text with your finger or stylus

8

To access other menus



Or



➔ Use this button to navigate between menus

9

Saving and loading images

Press



➔ To save image press



➔ To load an image press



10

Print

Press



➔ Sends image to printer

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
UGIC0709