


Upgrading the Firmware in a 4MPixel Chemi XT Camera

To upgrade the firmware in a GBox Chemi XT system:

1. Transfer the program **Atik_Firmware Uploader_4021_4_4.exe** to the computer connected to the camera.
2. Ensure that GeneSnap is NOT running.
3. Power on the darkroom and camera.
4. Run the program **Atik_Firmware Uploader_4021_4_4.exe**
5. When the introduction dialog appears press “Next” to continue.
6. When the dialog below appears it will show the current firmware version and the version of firmware that will be uploaded when you press the upload button. If the current firmware version is 4.3 or lower then press “Upload” to continue. If a dialog appears warning that the camera you are connected to doesn’t match the expected type then press “No” and contact support@syngene.com

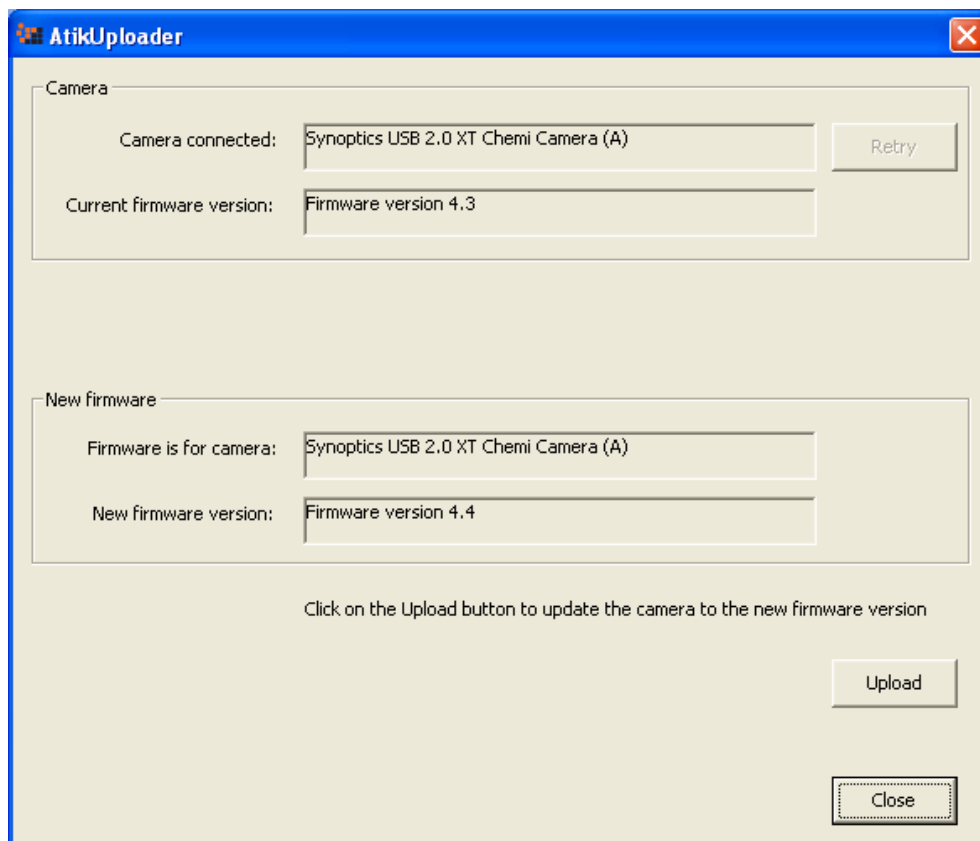

7. Uploading the new firmware will take just over 5 minutes. When complete the current version will show as version 4.4.
8. Press the close button to complete the upgrade.