

ChromaScan Image Capture-Quick Guide

1 Start ChromaScan

- ➔ Each user can enter a name or title for their own individual user preferences

2 Select configuration

- ➔ To set up system


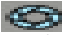

3 Live Image

Press




- ➔ The image will now be live

4 Position sample

- ➔ Open the cabinet and place the sample on the TLC viewing platform and close the door
- ➔ The user can adjust the Focus , Iris  and Zoom  settings if required.

5 Setting exposure time

- ➔ Click arrows to increase or decrease exposure time
- ➔ Or select Auto exposure box  This enables the software to automatically acquire a non-standard image

6 Grab image

Select



- ➔ The image will now be frozen

7 Save

Click



- ➔ Choose from Save As Or
- ➔ File/Export image to TIF, BMP, or JPEG

9 Print

Click



➔ Sends image to printer

10 Help menu

➔ Select the Help menu from the tool bar

➔ Use the drop down menu to pick a help topic

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

www.syngene.com
CSIC0709

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com