

DigiGenius Image Capture - Quick Guide

1 Turn on

Select 

➔ Turn on the camera (ensure the battery is charged)

➔ The Canon control software launches automatically when you turn on the camera. If the software does not start automatically click on Start/All Programmes/Canonutilities/EOS Utility/EOS Utility

2 Capturing an image

Select 

➔ Select Live View shoot to open live window

➔ Ensure the system has been installed correctly, as per the Installation Quick Guide

3 Settings

Select 

Select 

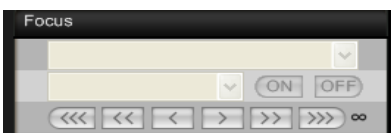
➔ From the White Balance menu select Cloudy from the drop down menu

4 Turn on transilluminator

Select 

➔ If bands on gel are faint then increase the white balance (From the EOS 1000D window right click on the control to activate it). For very faint bands go to +2

5 Focus setting

Select 

From the Remote Live View window

➔ To adjust zoom turn the lens of the camera

➔ To adjust focus use the arrows

6

Grab image

Press



From the EOS 1000D window

- ➔ The camera will set an exposure time relevant to the light levels coming from the sample
- ➔ Then turn off the camera

7

Save and export

- ➔ Images are saved as JPEG files
- ➔ Export as TIFF or BMP

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
DGIC0511