

GeneGnome⁵ Classic Image Capture - Quick Guide

1 Turn on

➔ Press the black button at the rear of the instrument

2 Start GeneSnap software

➔ Each user can enter a name or title for their own individual user preferences

3 Open drawer

Select



Open




Closed

➔ Place your sample in between the scored 8 position arrows present on the edges of the drawer frame. Then gently push the drawer to close

4 Choose appropriate settings

Select

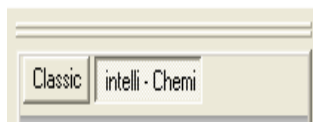

Press



Select




➔ Select the Classic tab
➔ When the door is shut the button will turn blue and the image is no longer live

➔ Press the blue button to start

➔ Select series capture to either capture a single frame or a series of images

➔ Using the drop-down menu select the number of images you wish to capture in the series

➔ Set Exposure time and click OK

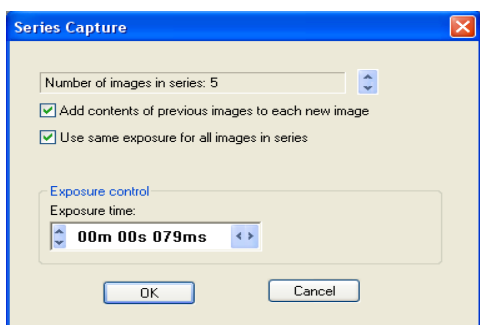

5

Save

Click



- ➔ Choose from Save As
- Or
- ➔ File/Export image to TIF, BMP, or JPEG

6

Print

Click



- ➔ Sends image to printer

7

Help menu

- ➔ Select the help menu from the toolbar
- ➔ Use the drop down menu to pick a help topic

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
GG5ClassicIC0211