

GeneGnome⁵ Installation - Quick Guide

■ System components

1. GeneGnome⁵ (red mini darkroom)
2. Integral PC
3. Flatscreen monitor with keyboard and mouse

■ Installation

Connecting the components

1. Unscrew the cover plate on the left-hand side of the GeneGnome⁵
2. Remove the cover plate to reveal the connectors beneath
3. Plug in the flat panel monitor, keyboard and mouse leads (the sockets are labelled with appropriate icons)
4. If a thermal printer has been supplied connect the printer cable to the appropriate socket
5. Refit the cover panel, slotting the cables through the cut-outs

CAUTION: Connect electrical power leads to each component but DO NOT SWITCH ON AT THIS TIME

■ Switching on

1. Ensure each mains lead is now connected to a live supply
2. Switch on the GeneGnome⁵ using the On/Off switch at the rear of the unit
3. Switch on the GeneGnome⁵ processor (hold the round button on the rear of the GeneGnome⁵ for 1-2 seconds). Also switch on the computer monitor and thermal printer (if connected)
4. The software is pre-loaded. You will need to accept the Windows licence

When the Windows operating system is fully loaded the GeneGnome⁵ system is ready to use

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
G65IC0211