

New Lighting Modules Installation Guide

■ Identifying modules

1. Each LED module is clearly labelled with a white sticker on the left hand side of the module.



■ Installation

Inserting modules

1. Remove any blank modules from LED holder.
2. Place LED module into LED module holder by placing a finger either side of the module.
3. Push the LED module in until you hear the module click into place.



LED
Module
Holder

Positioning modules

1. Modules of the same type i.e. two blue LEDs would normally be placed directly opposite each other.
- N.B. If you have a pair of LEDs make sure that you place the LED modules either side of the darkroom to ensure even sample illumination.

Removing modules

1. To remove a module place fingers top and bottom of the module pressing the clip on the bottom and gently glide out.

N.B. Be careful to not drop the module as this may damage it.

■ Transportation and storage of modules

1. The LED modules **MUST NOT** be inserted in the LED module holder during transit.
2. Wrap the LED modules in anti static bag or bubble wrap to prevent damage during transit and when modules are not in use.

CONTACT SYNGENE:

EUROPE:

Tel: +44 (0)1223 727123

Email: support@syngene.com

USA:

Tel: 800 686 4407/301 662 2863

Email: support@syngene.com

www.syngene.com

LMIG0710