

G:BOX Range Manual Series & Multiplex Image Capture - Quick Guide

1 Start GeneSys software

➔ Each user can enter a name or title for their own individual user preferences

2 Home screen

Select



➔ Select Manual Capture from the home screen

3 Position sample

➔ Place sample on UV transilluminator or visible light converter

N.B. If already in manual capture and the following controls (Lighting and filters) are not available then press the Home icon and then Manual capture

4 Capture Setup

Press



➔ Use Capture Setup to perform series capture and capture multiplexed images

Select



➔ Select to perform a Multiplex, Series capture or an additive series (contents of previous image added).

Select



➔ To perform a series capture select the Series button . Then select number of images you wish to capture then press the green arrow

Select



➔ If you wish to perform binning select the level of binning from the drop down menu

Select



➔ Select the appropriate lighting, filter and exposure time. If you wish to alter the lighting, filter or exposure time of one of the images in the series then press the green arrow then deselect same as previous

Uncheck

Same as previous

Press



➔ Press the green arrow to change between images then press capture

Press 

Select 

Select 

Select 

Select  

Uncheck Same as previous

Press 

➔ To perform a manual multiplex capture click Capture Setup

➔ Select Multiplex

➔ Select the number of images you wish to capture then press the green arrow

➔ If you wish to perform binning select the level of binning from the drop down menu

➔ Select the appropriate lighting, filter and exposure time. Then press the capture button

➔ Press the green arrow then deselect same as previous to select the appropriate lighting, filter and exposure time for each image. Press the green arrow to change between images

➔ Press the capture button

5 Print and save

Click 

➔ Click on the printer icon to send image to printer

Click 

➔ Click save and choose to save image as a TIF, BMP or JPEG.

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
GBRMSMI1110