


G:BOX Mini Monitor Mount Quick Guide

■ Components

1. 1 x G:BOX Mount bracket
2. 1 x Monitor Mount bracket
3. 1 x Monitor Mount spacer
4. 4 x M4x16 Screws
5. 8 x M4x12 Screws
6. 4 x M4x8 Screws

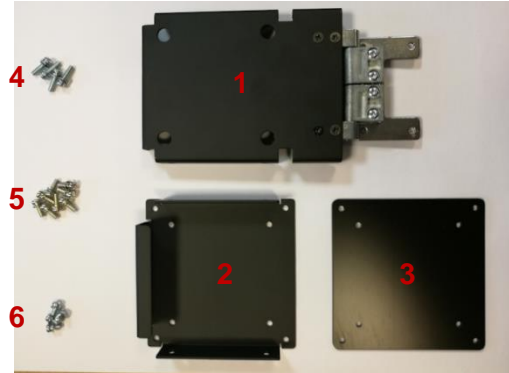

■ Attaching the mount

1. Remove the sticker from the front of the top cover.
2. Using 4 of the supplied M4x12 screws, attach the mount to the top cover as shown.
3. Carefully place your monitor face down on a clear surface taking care not to damage the screen. If the VESA mount on the monitor is recessed it is recommended that you use the supplied spacer between the monitor and the Monitor Mount bracket.
4. Using either the supplied M4x12 or M4x16 screws (dependant upon whether using the spacer or not), attach the Monitor mount (and possibly spacer) to the back of the monitor as shown.
5. Lower the monitor onto the G:BOX Mount bracket and secure at the sides of the bracket using the 4 M4x8 screws.



■ Additional Safety information

1. The G:BOX Mini mount is designed for up to 7kg. Any increase on this could cause damage to the top cover.

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
GBMM1117