

GeneSys StatusLink Broadcasting- Quick Guide

GeneSys can broadcast its status over a local network, allowing you to monitor the status of images you are capturing

1 Load GeneSys Software

- ➔ Each user can enter a name or title for their own individual user preferences

2 Set User Preferences

Select



- ➔ From the home screen select the 'Set User Preferences' icon
- ➔ Select the 'General Tab' on the left hand side of the screen

3 Network Broadcast Settings

Check

- ➔ In the 'Network Broadcast Settings' group check the 'Broadcast status over network' box
- ➔ The default port can be changed to do this. Please contact your network administrator to find out which ports you have access to

4 Save

Select



- ➔ Make sure that you press or select the 'Save' icon
- ➔ GeneSys will now register the requested port on your machine
- ➔ GeneSys will now display a pop-up message box giving the address of the page where it will broadcast its status over the network

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
GSSB1214