

NuGenius/NuGenius+ Installation Quick Guide

■ System components

1. Darkroom (Epi White lighting factory fitted)
2. High performance CCD camera with zoom lens pre-fitted to camera
3. UV filter (screwed on to the camera lens)
4. Full set of leads and cables

CAUTION: Do not connect power supply to any of the components until you are satisfied that everything is connected correctly. For assistance please contact your supplier or Syngene directly


■ Installation

Transilluminator set-up

1. Place the UV transilluminator inside and connect it to the free flying mains lead inside the NuGenius system
2. Turn on the transilluminator

Power on

1. Connect the mains lead from the NuGenius system to the mains power supply and switch on
2. After a few seconds a display will appear and the image capture software will automatically open
3. There is one LED at the side of the unit - the LED will remain on when the system is running to indicate that the system is on

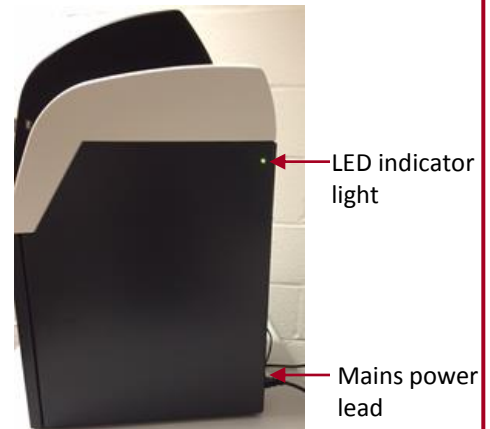

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
NGI1216