

U:Genius³ Installation - Quick Guide

■ System components

1. Darkroom (Epi White lighting factory fitted)
2. 3MP camera with zoom lens pre-fitted to camera
3. Stylus which can be attached to the darkroom in a suitable location
4. Full set of leads and cables

CAUTION: Do not connect power supply to any of the components until you are satisfied that everything is connected correctly. For assistance please contact your supplier or Syngene directly

■ Installation

Transilluminator set-up

1. Place the UV transilluminator inside and connect it to the free flying mains lead inside the U:Genius³ system

Camera set-up

1. Connect the USB cable from the camera to the back of the U:Genius³ unit

Installing the filter

1. Place the UV filter with the thread pointing upwards in the filter drawer (found below the camera and on the top of the U:Genius³ system)
2. This filter can be interchanged with other filters for different dyes



Power on/off

1. Connect the mains lead from the U:Genius³ system to the mains power supply and switch on
2. Power on **momentarily** pressing the round black button on the rear of the unit. There will be a 1-2 second delay as the system boots
3. After a few seconds a display will appear
4. To turn the internal computer off press the round black button on the rear of the unit



Select Regional Power Frequency

1. From the first screen select 'Preferences' icon
2. Select either 50 or 60Hz.



or



■ Installing GeneTools on a separate PC

1. Plug the Syngene branded flash drive into the PC you wish to install GeneTools on. Navigate into the "GeneTools" folder on the flash drive and run the "InstallGeneTools" program
2. Follow the on-screen instructions to install and license the GeneTools software. You will probably have been provided with media keys to license the software. These media keys can be found in a file named "Media Keys.txt" on the Syngene branded flash drive

CONTACT SYNGENE:
EUROPE:
Tel: +44 (0)1223 727123
Email: support@syngene.com

USA:
Tel: 800 686 4407/301 662 2863
Email: support@syngene.com

www.syngene.com
UG3INST0416